





Process Engineering
Division

Niro Pharma Systems

AEROMATIC
BUCK
COLLETTE
COURTOY
FIELDER
LYOPHIL
NICA
NIRO

Investigation of a New Coating Process for the Application of Enteric Coatings to Small Tablet Samples

A Summary of the
Poster Presented at the November
2005 AAPS Poster Session in
Nashville, TN

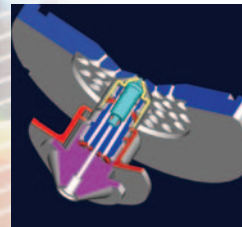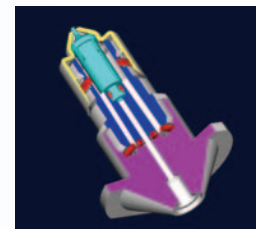

Investigation of a New Coating Process for the of Enteric Coatings to Small Tablet Samples

Charles Cunningham¹, Andrew Birkmire², Chris Neely¹ and Adam Kevra¹

Colorcon ¹Pharmaceutical Technical Services, West Point, PA, USA, GEA Niro Pharma System

Objectives The purpose of this study was to investigate the rapid application of an aqueous enteric coating system onto very small batch sizes of tablets using a novel coating process technology. The SUPERCELL™ (GEA Niro Pharma Systems) process involves rapid rotation of tablets as they pass continuously through the coating zone potentially allowing for very uniform film formation on all surfaces of the tablets. This technology could be advantageous where the quantity of active tablets is limited and not sufficient for successful coating in a rotating pan.

Experimental Methods



Aspirin tablets (325mg) were used as the coating substrate. An enteric coating system, Acryl-EZE® (Colorcon) was chosen for this evaluation since coated tablet defects can be quantitatively assessed using acid resistance testing. Even small defects in the tablet coating can result in failure of the enteric system.¹

Acryl-EZE is a fully formulated, pigmented aqueous enteric (methacrylic acid copolymer type C) coating system. The dispersion was prepared at 20% solids concentration. No subcoat was applied to the tablets prior to enteric coating.

Supercell process

In the Supercell process individual batches of 15 to 100 grams of tablets or capsules can be coated co-currently with heated drying air. The tablets are rapidly rotated as they pass through the coating zone ensuring intimate contact with the coating material (Figure 1).

Tablets are loaded into the chamber and conveyed through the coating zone in a ballistic flight path. Coating material is atomized by means of a specially designed low-momentum two-fluid spray nozzle located below the air distribution plate (Figure 2).

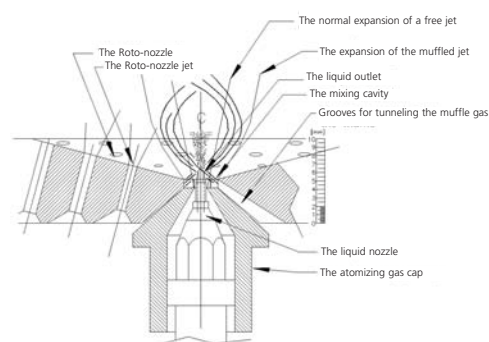


Figure 2: Schematic of spray nozzle and air distribution plate

The atomizing air is mixed with low-pressure drying air to dissipate the momentum generated in the atomizing process and reduce the potential for edge attrition of the tablets when striking the coater surfaces.² After coating, the tablets or capsules are pneumatically conveyed out of the coating chamber.

Six coating trials were conducted with batch sizes from 45g to 48g and target weight gains from 6 to 11% w/w. The coating machine automatically pre-weighed the tablets and fed them into the coating chamber with automated discharge at the completion of each coating cycle in a semi-continuous process. Batch recipes were programmed into the control system in advance of coating (Table1).



Figure 1: Tablet movement

References:

1. G.R.B. Down et.al., Drug Dev. Ind. Pharm. (19) page 2746, 1993
2. Birkmire et al., Poster presentation, AAPS National Meeting, Baltimore, 2004

Application

ms²Engineering Department, Columbia, MD, USA

| Trial | Airflow (m ³ /hr) | Inlet air temp. (°C) | Atomizing air pressure (bar) | Liquid volume applied (ml) | Spray rate (ml/min) |
|-------|------------------------------|----------------------|------------------------------|----------------------------|---------------------|
| 1 | 20.03 | 70 | 2.5 | 13.5 | 3.8 |
| 2 | 20.96 | 70 | 2.5 | 15.8 | 3.8 |
| 3 | 20.37 | 70 | 2.5 | 18.0 | 3.8 |
| 4 | 20.36 | 70 | 2.5 | 20.3 | 3.8 |
| 5 | 20.48 | 70 | 2.5 | 22.5 | 3.8 |
| 6 | 20.57 | 70 | 2.5 | 24.8 | 3.8 |

The resultant tablet samples were tested in 0.1N HCl for 2 hours for enteric properties using a modified disintegration apparatus.

In this modified disintegration method, 50 enteric-coated tablets from each trial were placed in a modified disintegration basket capable of holding 25-200 tablets. The tablets were subjected to 0.1N hydrochloric acid (37.5°C) for 2 hours. Results were reported as the percentage of tablets which exhibited no signs of cracking, peeling, bloating or disintegration after 2 hours.

Results

Tablet appearance

The coating process for each of the six trials ran smoothly with no difficulties. Visual examination of the tablets after coating revealed no signs of edge defects or other tablet damage. Edge defects were a concern due to random tablet impact with other tablets and various machine surfaces. The Acryl-EZE coated tablets exhibited exceptionally smooth and glossy surfaces (Figure 3).

Coating time

The tablets were exposed to the coating process for a very short period of time. The longest trial with a coating target of



Figure 3: Coated tablet appearance (10% weight gain)

11% weight gain was completed in 6.5 minutes. Trial 1 with a target weight gain of 6% was completed in 3.5 minutes. This short coating duration may have the potential to minimize the risks associated with temperature or moisture sensitive actives exposed to coating conditions for long periods of time.

Enteric testing

The enteric test results were very good with the tablets coated to a 9, 10, or 11% theoretical weight gain (trials 4-6) passing the enteric test requirements with less than 10% of the tablets exhibiting signs of defects. Trials 5 and 6 remained 100% intact after 2 hours exposure to acid (Figure 4). Tablet samples from trial 6 were tested for dissolution performance and met the USP criteria for delayed release aspirin.

Conclusions

The results indicate that the Supercell process enables a fast, uniform application of functional enteric coatings to very small batch sizes. The Acryl-EZE enteric coating system used in the study resulted in good enteric performance at application levels that would be typical in rotating pan technology. In early stage drug development where active materials may be in short supply, this coating technology will allow for evaluation of coating systems on a very small scale. It may also be possible to coat friable tablets successfully due to the fast development of the coating but care must be taken if scaling up the process to a rotating type coating pan. With the automated, semi-continuous operation of the machine and the very rapid coating application, it may be possible to link multiple machines together for full scale production of coated tablets.

Table 1: Main process parameters

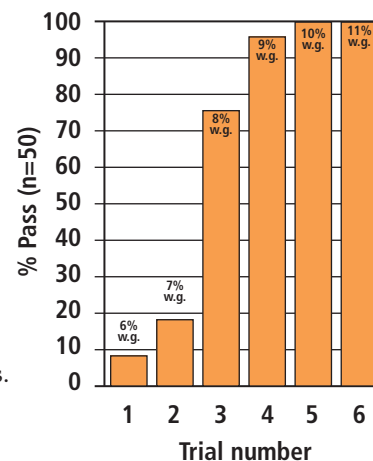


Figure 4: Enteric test results

Colorcon, Inc. is a global company that develops, manufactures and supplies specialty chemicals and excipients for the pharmaceutical industry. The core business is the design and technical support of advanced coating systems for pharmaceutical dosage forms for immediate, sustained and enteric release. In addition, the company's technological capabilities include assistance in the development of solid dosage formulations.

AEROMATIC
BUCK
COLLETTE
COURTOY
FIELDER
LYOPHIL
NICA
NIRO

Niro Pharma Systems is world leader in providing advanced processing solutions for solid dosage forms to the pharmaceutical industry. Based on a dedication to research and durable quality, Niro Pharma Systems offers a wide range of solutions, from individual pieces of equipment to complete integrated plants, by uniting the state-of-the-art technologies of Aeromatic, Buck, Collette, Courtoy, Fielder, Lyophil, Nica and Niro.

WWW.NIROINC.COM / WWW.COLORCON.COM



Niro Inc. • 9165 Rumsey Road, Columbia, MD 21045, USA • TEL +1 410 997 8700 • FAX +1 410 997 5021
E-MAIL info@niroinc.com • WEB www.niroinc.com

Colorcon, Inc. • 415 Moyer Blvd., P.O. Box 24, West Point, PA 19486-0024, USA • TEL +1 215 699 7733 • FAX +1 215 661 2605
E-MAIL info@colorcon.com • WEB www.colorcon.com